Model 1600

Neuroprobe Intracellular Amplifier





The Model 1600 Neuroprobe Amplifier is designed to meet the exacting requirements of intracellular recording and stimulation. The instrument consists of a high inputimpedance amplifier combined with current injection and bridge balance circuitry, allowing simultaneous stimulation and recording through the same electrode. The Model 1600 makes possible extremely precise recordings during current injection without requiring an external bridge or a differential input oscilloscope. A digital meter displays membrane potential, injected current, or electrode impedance.

The Neuroprobe Amplifier's advanced current injection system allows external switching of both internal and external current sources. The total current from all sources can be displayed on the digital meter prior to injection. An lontophoresis Adapter is available to allow for the application of high voltages for iontophoretic injection of drugs or dyes, or for any other application where greater current levels are required.

The Model 1600 Head Stage can be conveniently adapted to a wide variety of experimental situations and comes complete with a mounting rod and bracket for a standard micromanipulator and a standard 2 mm connector for a half-cell microelectrode holder.

Common applications for the Model 1600 Neuroprobe Intracellular Amplifier include, but are not limited to:

- o Intracellular recording of single cell action potentials
- o Current Clamp
- Dye Injection
- o Ultra-low noise extracellular recordings

The Model 1600 Neuroprobe Intracellular Amplifier has been an industry leader for over 25 years, and is still among the quietest recording amplifiers available on the market.



x10 Headstage Probe

- Digital display of membrane potential, injected current, or electrode resistance
- Square-wave generator for testing electrode resistance and capacitance compensation
- Current injection system adjustable to 0.1 nA resolution
- Simultaneous stimulating and recording by a single electrode
- Transient and DC balance controls
- Low-pass and Notch filters
- Includes rack mount hardware
- 3-year warranty

Model 1600

Neuroprobe Intracellular Amplifier



Application **Notes:**

extracellular recording of action potentials

While the Model 1600 Neuroprobe Intracellular Amplifier was designed to record intracellularly, many researchers utilize the Model 1600 in their extracellular recordings. The high input impedance of the amplifier results in extremely low noise recordings. The output of the Model 1600 Neuroprobe Amplifier is then routed to an additional gain amplification stage and perhaps a high pass filter before ultimately being recorded by a digitizing data acquisition system. For more information, contact A-M Systems or your distributor.

Specifications

10¹¹ ohms Current Input Impedance

Current Input Capacitance Adjustable to zero

Current Input Range +/-10 V; Up to 200 V when using the optional lontophoresis Adapter

Current Injection 10 nA / V or 100nA / V Input Bias Current Adjustable to zero Frequency Response DC to 325 kHz

Low Pass Filter 1, 2, 5, 10, 20, 50, and 100 kHz; -12 dB / octave

Notch Filter 50 Hz or 60 Hz Capacity Compensation -4 pF to 30 pF

Noise 0 ohm source: 13 microvolts rms (10 Hz to 50 kHz)

> 1 megohm source: 79 microvolts rms (10 Hz to 50 kHz) 20 megohm source: 310 microvolts rms (10 Hz to 50 kHz)

DC Balance Low range: up to 500 megohms

High range: up to 50 megohms

Outputs x1 or x10 gain

Current Monitor Low range: 10 mV / nA

High range: 1 mV / nA

References

Mason AC, Oshinsky ML, and Hoy, RR (2001) Hyperacute directional hearing in a microscale auditory system. Nature 410: 686-690

Person AL and Perkel DJ (2007) Pallidal neuron activity increases during sensory relay through thalamus in a songbird circuit essential for learning. J Neuroscience 27(32):8687-8698

Yoon AC et al. (2008) Determining CA^{2+} -sensor binding time and its variability in evoked neurotransmitter release.

Ordering Information

Product #680105 Country-specific power cords are not supplied. For use on 220 V / 50 Hz power systems:

For use on 110 V / 60 Hz power systems: Product #680100 Product #682000 Optional: Iontophoresis Adapter:

One product #681500 Headstage probe must be purchased at time of ordering for proper operation of the

instrument.

J Physiology 586:1005-1015

All units include a product manual and rack mounts.

Distributed By:

Science Products GmbH

Hofheimer Straße 63 D- 65719 Hofheim

Phone: +49 6192 90 13 96 Fax: +49 6192 90 13 98

Science Products Trading SPT AG

Käferholzstr. 142 CH- 4058 Basel

in collaboration with:

Phone +41 43 488 05 61 Fax: +41 43 488 05 62

Email: info@science-products.com

+48 12 363 80 52

+43 7207 7 54 55